Surface Concept Devices

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Safety Instructions (GB, USA)

Surface Concept devices operate with lethal high-voltages.

All Surface Concept safety instructions must be strictly followed.

Surface Concept safety instructions should always be kept near the device, so that they can be easily and quickly accessible for all persons.

Correct operation and use of the device only for its intended purpose will ensure that the device will give trouble-free performance and will remain in safe condition.

Regulations for protection against radiation sources must also be observed.

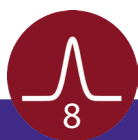
Important Information

For more detailed instructions on maintenance and for the requirements governing safety of the device, the relevant national, resp. federal regulations are to be noted and strictly observed.

If these regulations are not known, a specialist should be consulted in all cases!

Surrounding Conditions

- After the device is unpacked, it must be allowed to sit for at least 24 hours at the installation location (possible formation of condensation). Only then should the power connection be made.
- The device must be operated in a dry, covered area.
- The device must be placed on a flat, level and no-slip work surface, and be fitted into a suitable 19" rack, where appropriate.
- The device must not be exposed to extremely high or low temperatures.
- The device must not be exposed to direct sunlight and not be stored, installed or operated close to heating systems, heat generating devices and/or air-conditioners.

- The permissible range for operation temperature is defined between 5°C and 40°C.
- The device must not be operated in areas of high humidity (for example frost, dew, condensation or rain water), draught, dust accumulation or danger from oil splashes, iron dust, and corrosive liquids.
- The maximum relative air humidity is defined at 80% at 31°C (50% at 40°C).
- The device must be placed so that the air ventilation is not obstructed. The specific guidelines in the operation instructions are to be followed.
- Trouble-free operation can be guaranteed for an altitude of up to 2000 meters above sea level.

Danger for Pacemakers

Persons with cardiac pacemakers and other surgical implants should not expose themselves to electromagnetic fields unless their pacemakers or implants comply with DIN VDE 0750 section 9A, EN 50061, ISO 601 -2-31 * CEI.

Main Connections

- Care is to be taken that the mains voltage agrees with the specifications on the identification plate on the rear of the device.
- Mains power supply-voltage fluctuations must not exceed +/- 10% from the rated value.
- The device must under no circumstances be connected with mains power supply outlets already used by other devices which could (possibly) cause interference.
- Before the device is switched on, all plugs must be securely inserted into their corresponding jacks. Screw connections are to be securely tightened and bayonet connections must lock in.
- If the device emit sparks, smokes, or smells of burning, it must be switched off immediately as well as all peripheral devices attached.
- In such an event any EMERGENCY OFF device fitted should also be immediately engaged.
- **Never use water to extinguish fire!**

Operation

- Before the device is started, the operator must ensure that it is in the proper condition and operating safely.
- Do not remove safety fittings, such as covers, screw joints, locking wires and seals for/before operation.
- Use the device only for its intended purpose.

Switching Off

- After the device has been switched off, the capacitors may continue to be charged with lethal voltages for some time. Check the voltage in all capacitors before any maintenance work is commenced.
- Work with radiation sources, cables, connected peripheral devices or cables only when the sources and all peripheral devices have been switched off.
- After the device has been switched off, cable connections may be removed only after a waiting period of one minute has elapsed.

Opening A Device

- The device may only be opened by specialists with detailed knowledge of the devices/ systems on the basis of special training and experience.
- Before the device is opened, all terminals must be disconnected from the power supply: Press the power switch (OFF or O Position) and remove the power plug (only on plug casing).
- In the case of devices with voltage over 5 kV, covers are only to be opened with suitable tools. Care is to be taken that only non-conductive, high voltage tools must be used.

Never put a device into operation when the casing is open!

Maintenance

- Devices, connections, power plugs and measuring equipment are to be subject to inspection for noticeable defects every six months.
- Cover all high-voltage conducting parts during maintenance work on opened devices.
- Acoustic and visual warning signals in the area of the devices/ units/systems should be observed and acted upon.
- As protection against hazardous electric shocks, external earthing equipment (twisted copper wire of 10mm to 15mm in width) must be subject to regular mechanical and electrical inspections for corrosion.
- The ground wire should be examined for power resistance every three months.
- **If the device/system is operating in an unknown operating condition, it should only be approached with the utmost caution, e.g.: With the left hand in the trouser pocket or behind the belt, the right hand carefully approaches the device.**

Cleaning

- The device may only be cleaned by qualified personnel.
- Before cleaning the device, care is to be taken that all peripheral devices are switched off and all plugs are removed from the jacks (complete disconnection of mains terminals).
- Only a soft dust-free cloth may be used. The cloth should be lightly moistured with water or a mild, non-aggressive liquid detergent (observe instructions for use).
- Specific care is to be taken that no moisture enters the inside of the device.

Never use strong chemical agents, such as alcohol, paraffin or paint thinner.

Accessories

- Only use original cables, plugs and accessories manufactured by Surface Concept.
- These components are to be subject to regular tests. Care is to be taken that they show no evidence of any mechanical and/ or electrical defects.
- In case of doubt, they are to be replaced immediately by original Surface Concept components.
- When fuses are being replaced, only the same type must be used: voltage strength, strength of rated current and switch characteristics.

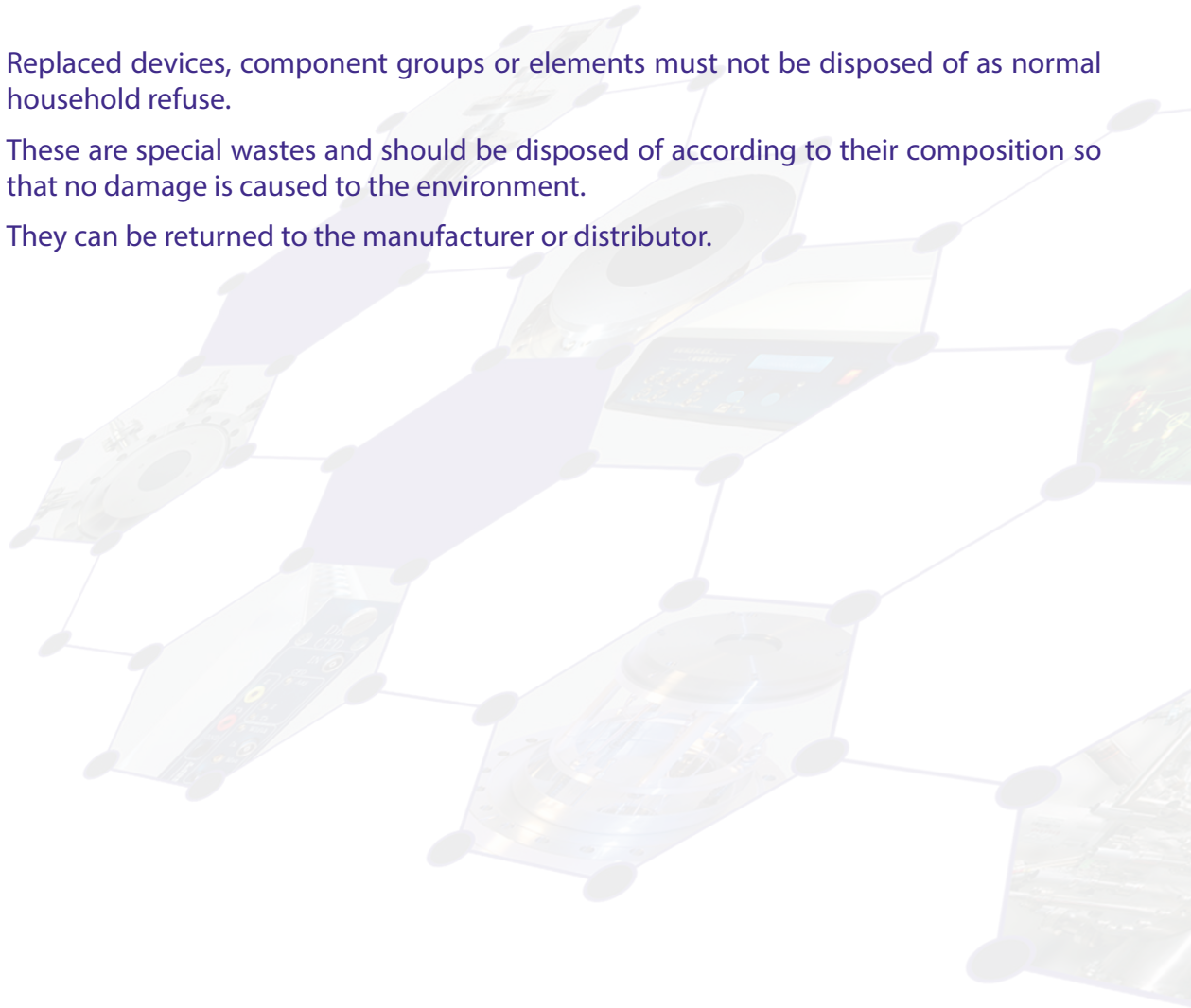
- If cables or other devices not manufactured by Surface Concept are connected and put into operation or used, there can be no warranty for the correct function for connected devices and systems nor for the safety of personnel.
- If there is any doubt, Surface Concept should be consulted immediately.

Protection of the Environment

- Replaced devices, component groups or elements must not be disposed of as normal household refuse.
- These are special wastes and should be disposed of according to their composition so that no damage is caused to the environment.
- They can be returned to the manufacturer or distributor.

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All rights reserved. No part of this manual may be reproduced without the prior permission of Surface Concept GmbH.